AquaCHROM™

Chromogenic medium for detection of *E.coli* and coliforms in water samples.

article number AQ056 : 100 x 100 ml doses

STORAGE :

Store the product at 15/30°C until the shelflife date indicated on the label.

COMPOSITION in g/L :

Peptones and growth regulators 20,0 ; Chromogenic and selective mix 2,3 ; pH : 7,1 +/- 0,2 (Classical formula adjusted and/or supplemented as required to meet performance criteria).

INOCULATION and INTERPRETATION :

- 1 Add a pre-weighed dose of medium in a sterile transparent vessel containing the 100ml water sample.
- 2 Close and shake until dissolution.
- 3 Incubate at 35-37°C for 18-24h.
- 4 Read the results (Presence/Absence) :

Typical appearance after incubation	Reading
-------------------------------------	---------

Green to blue-green	→ presence of <i>E.coli</i>
Yellow	→ presence of non- <i>E.coli</i> coliforms
Colourless ± growth *	→ other

* This selective medium is inhibitory for many microorganisms, mostly gram positive.

PERFORMANCE and LIMITATIONS :

- Some rare strains of *E.coli*, like *E.coli* O157, may display yellow colouration.
- Some rare strains of *Salmonella* can give a dark blue colouration.
- Colourless results can be obtained with some atypical strains of coliforms.

DISPOSAL OF WASTE :

After interpretation all media should be destroyed by autoclaving at 121°C for at least 20 minutes.

English. For laboratory use. Laboratory product to be used only by trained personnel.

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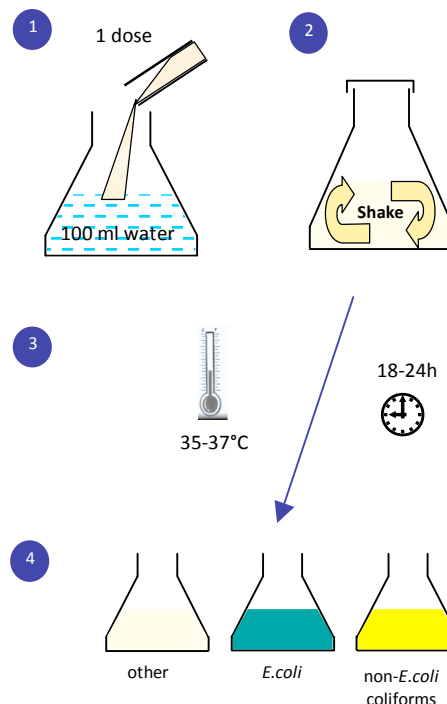
AquaCHROM Procedure

Available from CHROMagar :

CHROMagar™ Candida
Differentiation of major pathogenic *Candida* species

CHROMagar™ Orientation
Differentiation of urinary tract pathogens

Rambach™ Agar
Detection of *Salmonella* spp

CHROMagar™ Salmonella
Detection of *Salmonella* including *S. Typhi*

CHROMagar™ Salmonella Plus
Detection of *Salmonella* according to the ISO 6579:2002 norm

CHROMagar™ O157
Detection of *E.coli* O157

CHROMagar™ E.coli
Detection and enumeration of *E.coli*

CHROMagar™ ECC
Detection and enumeration of *E.coli* and coliforms

CHROMagar™ Liquid ECC
Broth for pad technique for *E.coli*-coliforms

CHROMagar™ Staph aureus
Detection and enumeration of *Staphylococcus aureus*

CHROMagar™ MRSA
Detection of MRSA including low level MRSA

CHROMagar™ Listeria
Detection and enumeration of *Listeria monocytogenes*

CHROMagar™ Vibrio
Detection and enumeration of *Vibrio parahaemolyticus*, *Vibrio vulnificus* and *Vibrio cholerae*

CHROMagar™ VRE
Detection of *E.faecium* VRE & *E.faecalis* VRE

CHROMagar™ StrepB
Detection of *Streptococcus agalactiae* (GBS)

CHROMagar™ ESBL
Detection of ESBL-producing bacteria

CHROMagar™ KPC
Detection of Carbapenem-resistant bacteria

CHROMagar™, Rambach™ and AquaCHROM™ are trademarks of Dr. A. Rambach

Visit CHROMagar on internet via <http://www.chromagar.com>

CHROMagar
Microbiology

4, place du 18 Juin 1940
75006 Paris France
Fax: (33-1) 45 48 06 06

NT-EXT-036	Version 2 30-Dec-2010
------------	--------------------------