CHROMagar™ B.cereus

Chromogenic medium for the detection and enumeration of *Bacillus cereus* group

Composed by:
CHROMagar™ B.cereus base: article number BC732(B)
CHROMagar™ B.cereus supplement: article number BC732(S)
(Sold as a pack under article number BC732)

STORAGE

Store the powder base at 15/30°C and the supplement at 2/8°C until the shelflife date indicated on the label.

COMPOSITION in g/L

Base: Agar 15.0; Peptone and yeast extract 8.0; NaCl 10.0; Chromogenic mix 0.3.

Supplement: 3.0.

→ pH : 6,8 +/- 0,2 (Classical formula adjusted and/or supplemented as required to meet performance criteria).

PREPARATION (for 1L of final media) :

• BASE :

Suspend 33.3 g of the CHROMagar B.cereus base into 1L of purified water. Disperse powder slowly in water by rotating for swelling of the agar. Heat to 121°C +/- 1°C during 15 minutes. Cool in a water bath to 47°C +/- 2°C.

• SUPPLEMENT :

Disperse slowly 3g of CHROMagar B.cereus supplement in 40ml of purified water. Add a magnetic bar and homogenize during at least 30 minutes rotating the magnetic bar at high speed (~1200 rpm) until obtaining a creamy homogeneous suspension. Heat to 121°C +/- 1°C during 15 minutes. Cool in a water bath to 47°C +/- 2°C.

• MIXING OF BASE AND SUPPLEMENT :

Place the melted 47°C cooled base under gentle stirring. Add aseptically the 40ml of the homogeneous reconstituted supplement, keeping the gentle stirring during 1 or 2 minutes until complete homogenisation. Pour IMMEDIATELY into sterile Petri dishes and let them cool. Let dry. Store in the dark.

Prepared plates can be kept for one day at ambient temperature, or for up to two weeks at (2/8°C) if properly prepared and protected from light and dehydration.

INOCULATION :

If the agar plate has been refrigerated, allow to warm to room temperature before inoculation. Streak or spread sample onto plate and incubate at 30°C for 18-24 hours.

INTERPRETATION

Microorganism → Typical colony appearance

***Bacillus of cereus* group** → blue, with white halo

Other *Bacillus* → blue, colourless, or inhibited

Gram negative bacteria → inhibited

Yeast and moulds → inhibited

LIMITATIONS

• Depending of the strains, morphology of the colonies of *Bacillus cereus* group could vary. As an example, *B.mycoides* typically displays rhizoid form colonies.
• Definite identification may require additional testing, as those described in the ISO 7932 norm.

DISPOSAL OF WASTE

After interpretation all plates should be destroyed by autoclaving at 121°C for at least 20 minutes.

English. For laboratory use. Laboratory product to be used only by trained personnel.

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CHROMagar™ and Rambach™ are Trademarks of Dr. A. Rambach