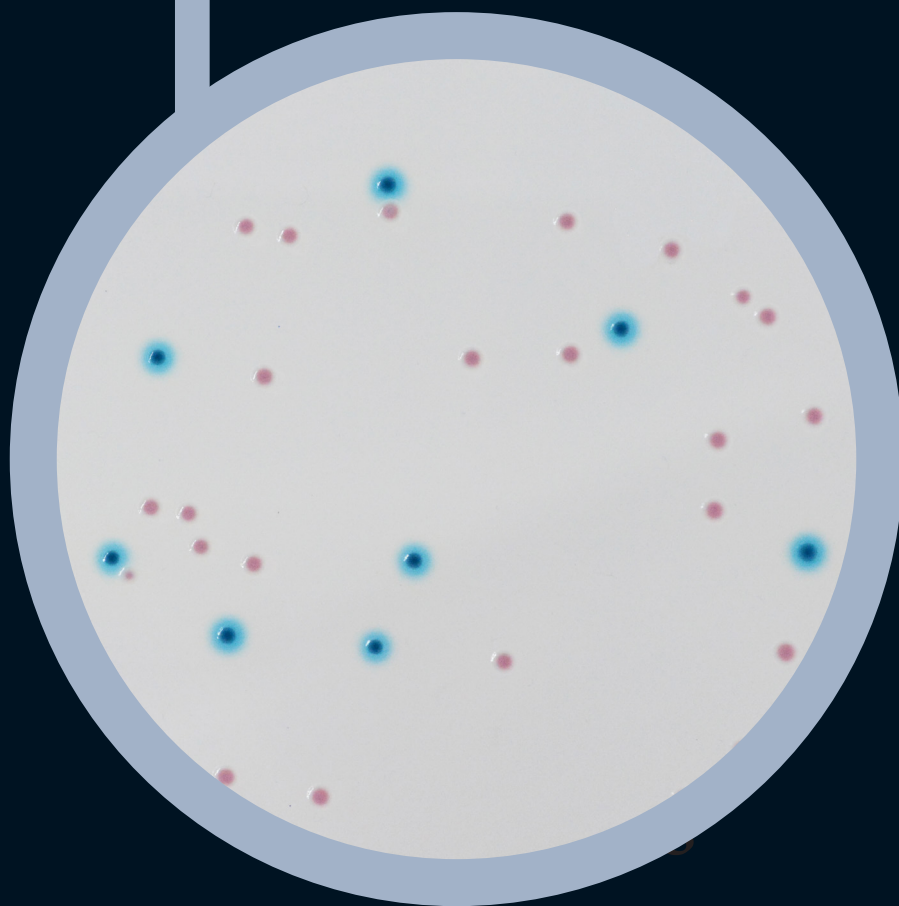
● CHROMagar™ Enterobacteria

**For detection and enumeration of
*Enterobacteriaceae***

● CHROMagar™ Enterobacteria


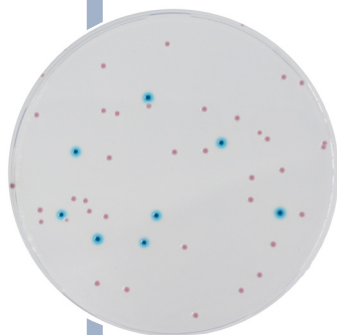
Plate Reading

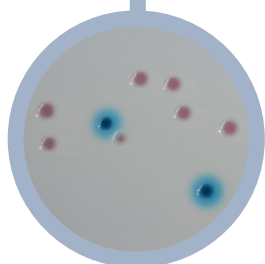- *E. coli*
→ blue with or without blue halo
- Other *Enterobacteriaceae*
→ pink to red
- Gram positive
→ inhibited

For detection and enumeration of *Enterobacteriaceae*

Background

«The *Enterobacteriaceae* and coliform bacteria within this family represent two of the most common groups of indicator organism used by the food industry. In some countries, depending on regulatory requirements, the food industry has moved towards testing for *Enterobacteriaceae*.» ILSI Europe (International Life Sciences Institute)

The ISO 21528 specifies a method for the enumeration of *Enterobacteriaceae*, applicable to products intended for human consumption, feeding of animals, and environmental samples using VRBG as culture medium.

CHROMagar™ Enterobacteria allows the detection and differentiation by the color of *E. coli* and other *Enterobacteria*.

Medium Performance

1 **EASY DIFFERENTIATION BETWEEN *E. COLI* AND OTHER *ENTEROBACTERIACEAE***
Unlike with VRBG, with CHROMagar™ Enterobacteria you can easily differentiate *E. coli* from the rest of *Enterobacteriaceae* by the color.

2 **EASY READING**
The intense colored colonies on a clear agar background facilitate the reading, contrary to VRBG where the smoked background allows little contrast.

Medium Description

Powder Base	Total	42.3 g/L
	Agar	11.0
Optional Supplements for ready-to-use bottles (to be ordered separately)	Peptones and Yeast extract	22.0
	Mineral salts	4.9
	Chromogenic and selective mix	0.7
	Growth factors	3.7
	Storage at 15/30°C - pH: 7.4 +/- 0.2	
	Shelf Life	2 years
	1 st : Liquid form.....	1 ml/L
Storage at 2-8 °C	Shelf Life	2 years
Usual Samples	2 nd : Powder form qsf 500 mL	0.01 g/L
	Storage at 2-8 °C	Shelf Life
Procedure	Food, foodstuff, environmental samples	
	Pouring, streaking or spreading techniques. Incubation 24h, 30 °C or 37 °C. Aerobic conditions.	

Order References

Please use these product references when contacting your local distributor:

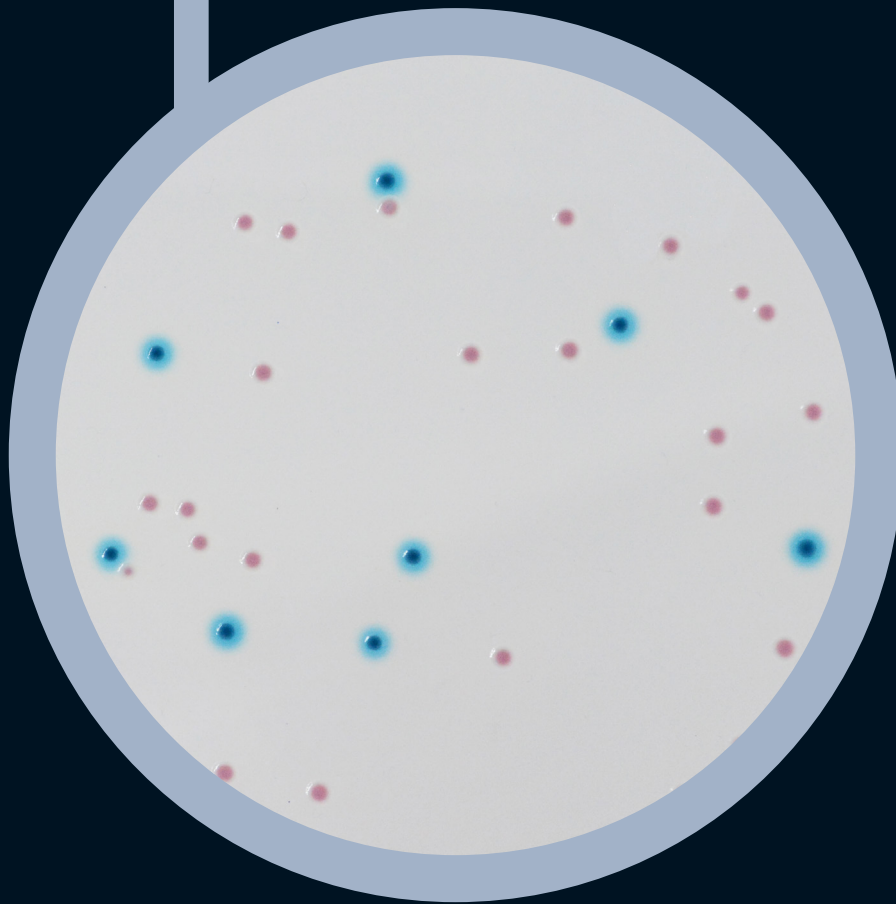
5000 ml pack EB042

Supplement 1 (to be used only with ready-to-use bottles)..... TT100L

Supplement 2 (to be used only with ready-to-use bottles)..... CV702

Manufacturer: CHROMagar
4 place du 18 juin 1940 75006 Paris - France
Email: CHROMagar@CHROMagar.com
Website: www.CHROMagar.com
Find your nearest distributor on
www.CHROMagar.com/contact

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