# AquaCHROM™ ECC

## Instructions For Use

Available in several languages

**NT-EXT-36**

Version **6.0**

Click below for:

**ENGLISH**

**FRANCAIS**

**ESPAÑOL**

## MEDIUM PURPOSE

Chromogenic medium for the detection of *E. coli* and coliforms in water samples.

Coliforms, *Enterobacteriaceae* able to ferment lactose, are bacteria present in human and warm blooded animals intestinal flora, in the soil and water. Coliforms are indicators of organic, environmental or faecal contamination.

Strict regulations exist for *E. coli*/coliform absence in water and food samples. This can be explained by the importance of these germs in determining water and food safety.

Worldwide, water and food quality control for human consumption are based on detecting the presence/absence of *E. coli* and coliforms.

## COMPOSITION

The product is composed of a powder medium.

Product	=	Pack
Total g/L		22.3 g/L
Composition g/L		Peptone and growth regulators 20.0 Chromogenic and selective mix 2.3
Aspect		Powder Form
<b>STORAGE</b>		<b>15-30 °C</b>
<b>FINAL MEDIA pH</b>		7.1 +/- 0.2

## PREPARATION

### Preparation

- This product is supplied ready to use, in pre-weighed doses.
- Each dose is for a 100 mL water sample.

### Storage

- Store in the dark before use.
- The product can be stored until the shelf life indicated on the label.

## INOCULATION

Water samples are to be processed as follows:

1. Add a pre-weighed dose in a sterile transparent vessel containing the 100 mL water sample

2. Close and shake until dissolution

18-24 h 35-37 °C

3. Incubate at 35-37 °C for 18-24 h.


4. Read the results (*E. coli*/coliforms).

## INTERPRETATION

Microorganism	Typical appearance
<i>E. coli</i>	→ green to blue-green
Other coliform bacteria	→ yellow
Other bacteria	→ colourless or inhibited

### Typical appearance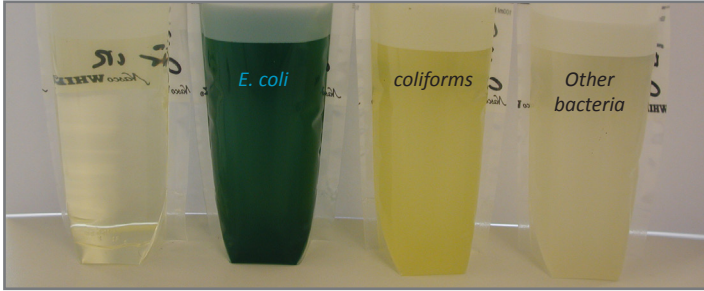

The pictures shown are not contractual.

## PERFORMANCE & LIMITATIONS

- Some rare strains of *E. coli*, like *E. coli* O157, may display yellow coloration.
- Some rare strains of *Salmonella* can give a dark blue colouration.
- Colourless results can be obtained with some atypical strains of coliforms.

## QUALITY CONTROL

Please perform Quality Control according to the use of the medium and the local QC regulations and norms. Good performance of the medium can be tested, inoculating the ATCC strains below:

Microorganism	Typical appearance
<i>E. coli</i> ATCC® 25922	→ green blue
<i>C. freundii</i> ATCC® 8090	→ yellow
<i>S. aureus</i> ATCC® 25923	→ inhibited

## WARNINGS

- Do not use the product beyond its expiry date or if it shows any evidence of deterioration.
- For Laboratory use. This laboratory product should be used only by trained personnel in compliance with good laboratory practices.
- Any change or modification in the procedure may affect the results.
- Any change or modification of the required storage temperature may affect the performance of the product.
- Unappropriate storage may affect the shelf life of the product.
- For a good microbial detection: collection and transport of specimen should be well handled and adapted to the particular specimen according to good laboratory practices.

## DISPOSAL OF WASTE

After use, all containers and any other contaminated materials must be sterilized or disposed of by appropriate internal procedures and in accordance with local legislations. Containers can be destroyed by autoclaving at 121 °C for at least 20 minutes.

## REFERENCES

Please refer to our website page «Publications» for scientific publications about this particular product.

Web link: <http://www.chromagar.com/publication.php>

## IFU/LABEL INDEX

- Catalogue reference
- Consult instructions for use
- Quantity of powder sufficient for X liters of media
- Expiry date
- Required storage temperature
- Store away from humidity
- Protect from light
- Manufacturer

Pack Size

100 doses of 100 mL

=

Ordering References

**AQ056**

Weight: 100 x 2.23 g

### Need some Technical Documents?

Available for download on [www.CHROMagar.com](http://www.CHROMagar.com)

- Certificate of Analysis (CoA) --> One per Lot
- Material Safety Data Sheet (MSDS)

CHROMagar™ and Rambach™ are trademarks created by Dr A. Rambach  
ATCC® is a registered trademark of the American Type Culture Collection  
NT-EXT-036 V6.0 / 26-Nov-18

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