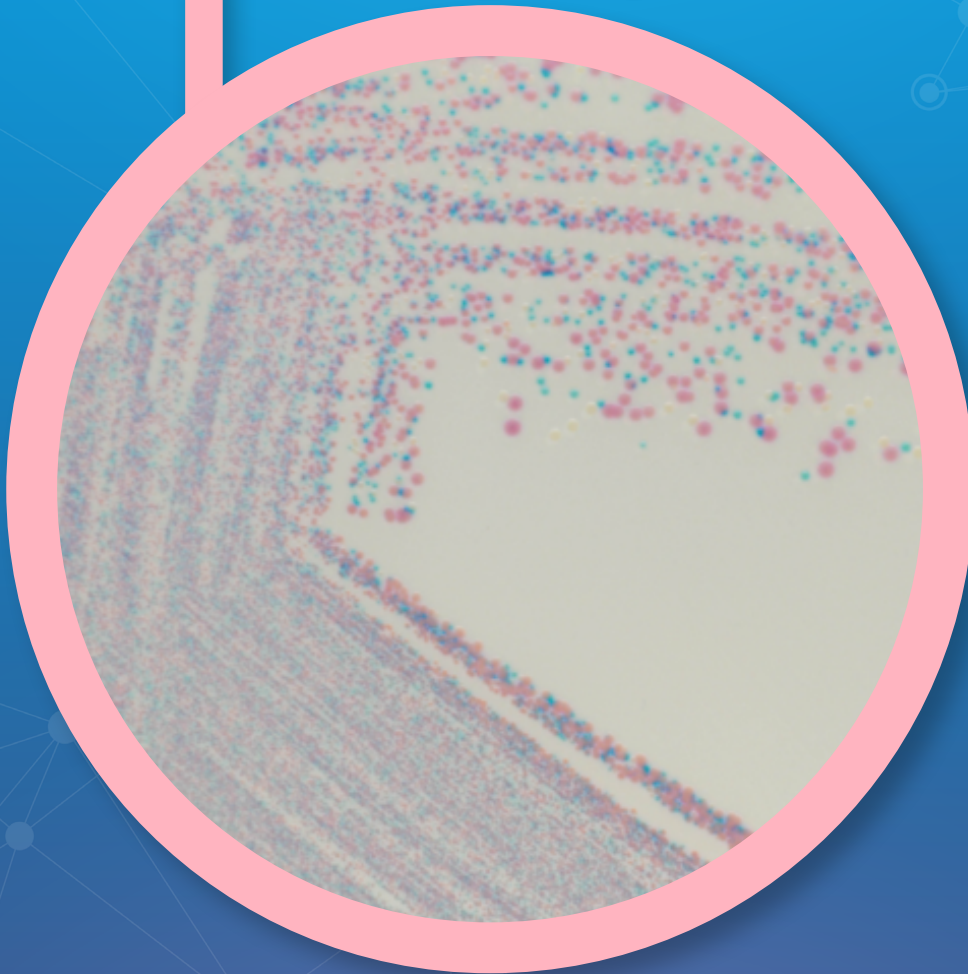
Leaflet

LF-EXT-049

Version 2.0

● CHROMagar™ Staphylococcus

Click below:

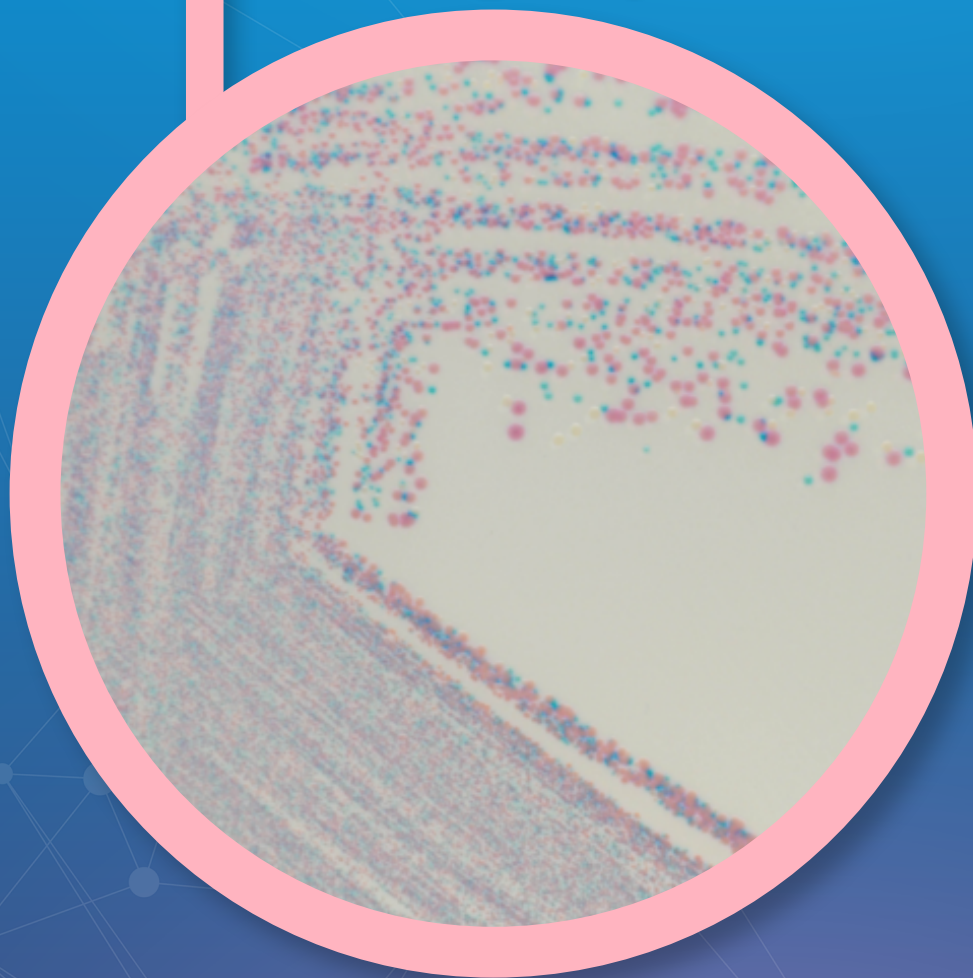
EN

FR

ES

CHROMagar™
The Chromogenic Media Pioneer

● CHROMagar™ Staphylococcus

**For isolation and differentiation
of *Staphylococcus* spp.**

● CHROMagar™ Staphylococcus

www.CHROMagar.com

For isolation and differentiation of *Staphylococcus* spp.


Background

Mastitis is a disease that causes major losses in dairy production. The most prevalent form of the disease is subclinical mastitis, which leads to decreased milk production and an increased risk of clinical mastitis during lactation. Controlling the spread of contagious subclinical mastitis by rapid identification of causing pathogens such as *Staphylococcus aureus* and other *Staphylococcus* species is fundamental for rapid control measures.

CHROMagar™ Staphylococcus is a selective and differential chromogenic medium for the qualitative direct detection and differentiation of *Staphylococcus* spp. In combination with CHROMagar™ Streptococcus and CHROMagar™ Mastitis GN, CHROMagar™ Staphylococcus becomes a powerful tool to aid in the detection of pathogens in milk samples. Results can be interpreted after 24 h of aerobic incubation at 37 °C.

Plate Reading

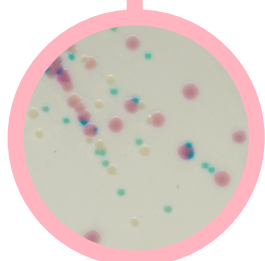- *S. aureus*
→ pink to mauve
- *S. epidermidis*
→ colourless to pinkish
- *S. saprophyticus*
→ turquoise blue

Medium Performance

1 EASY TO PREPARE

CHROMagar™ Staphylococcus comes with all compounds already included in the powder (no supplement needed), while the conventional medium for *Staphylococcus* spp., the Baird-Parker medium, must be supplemented with an emulsion of egg yolk and antibiotics after autoclaving, adding steps to plate manufacturing. CHROMagar™ Staphylococcus powder as well as the poured plates are stable for an extended shelf-life period.

2 FAST AND EASY READING

The results on CHROMagar™ Staphylococcus plates are available already after 24 h, while results are read after 48 h of incubation on Baird-Parker plates. In addition, CHROMagar™ Staphylococcus enables direct differentiation of *Staphylococcus* species by coloration and prevents growth of the unwanted annex flora through an enhanced selectivity.

Medium Description

Powder Base	Total 87.4 g/L Agar 15.0 Peptone and yeast extract 40.0 Salts 30.0 Chromogenic mix 2.4 Storage at 15/30 °C - pH: 7.3 +/- 0.2 Shelf Life > 18 months
Usual Samples	Milk, surfaces
Procedure	Direct streaking. Incubate at 37 °C for 24 h. Aerobic conditions.
Scientific Publications on this product: available on www.CHROMagar.com Please read carefully the instructions for use (IFU document) available on www.CHROMagar.com	

Order References

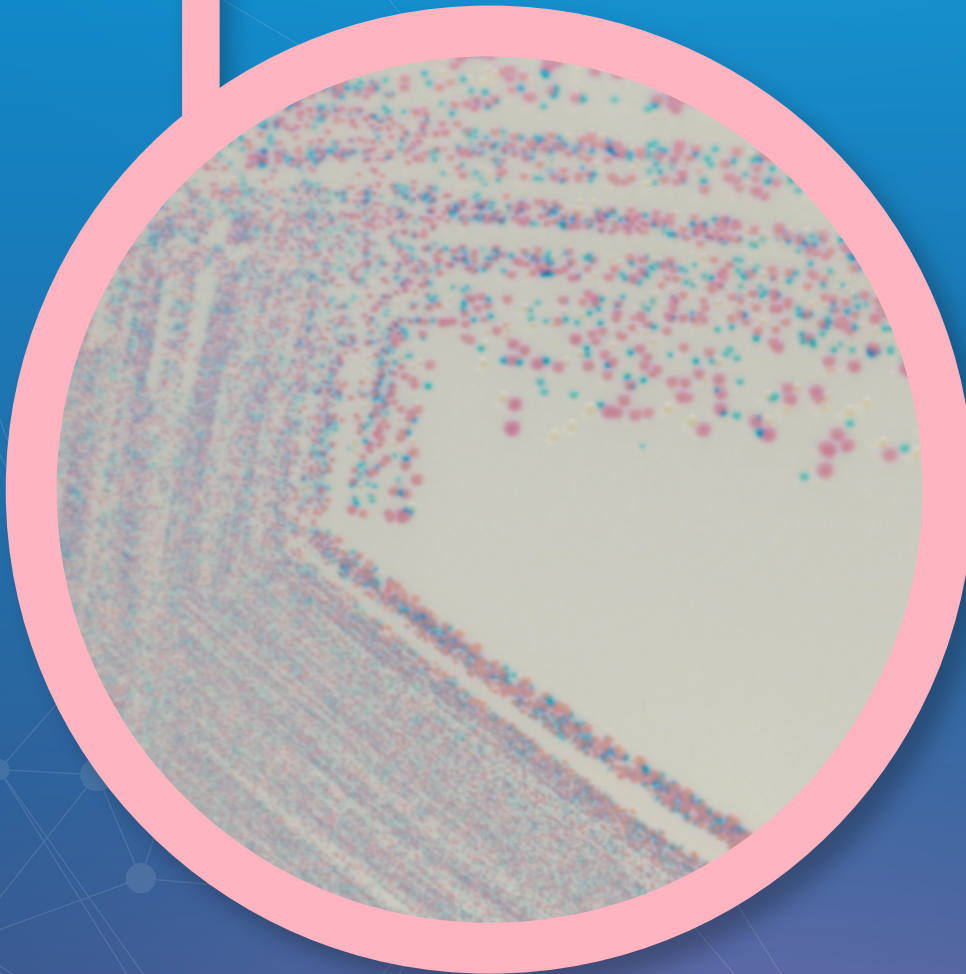
Please use these product references when contacting your local distributor:

500 g pack CQ38-500G

Manufacturer: CHROMagar, 29 avenue George Sand, 93210 La Plaine Saint-Denis - France
Email: CHROMagar@CHROMagar.com
Website: www.CHROMagar.com

Find your nearest distributor on
www.CHROMagar.com/contact

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