## CHROMagar™ Enterobacteria

Click below:

# CHROMagar™ Enterobacteriaceae

For detection and enumeration of  
*Enterobacteriaceae*


**CHRO** Magar™  
The Chromogenic Media Pioneer

# ● CHROMagar™ Enterobacteria

## Plate Reading

- *E. coli*  
→ blue with our without blue halo
- Other *Enterobacteriaceae*  
→ pink to red
- Gram (+)  
→ inhibited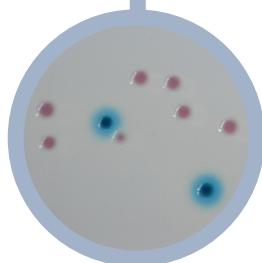

## For detection and enumeration of *Enterobacteriaceae*

### Background

«The *Enterobacteriaceae* and coliform bacteria within this family represent two of the most common groups of indicator organism used by the food industry. In some countries, depending on regulatory requirements, the food industry has moved towards testing for *Enterobacteriaceae*.» ILSI Europe (International Life Sciences Institute)

The ISO 21528 specifies a method for the enumeration of *Enterobacteriaceae*, applicable to products intended for human consumption, feeding of animals, and environmental samples using VRBG as culture medium.

CHROMagar™ Enterobacteria allows de detection and differentiation by the color of *E. coli* and other Enterobacteria.

### Medium Performance

#### 1 EASY DIFFERENTIATION BETWEEN *E. COLI* AND OTHER *ENTEROBACTERIACEAE*

Unlike with VRBG, with CHROMagar™ Enterobacteria you can easily differentiate *E. coli* from the rest of *Enterobacteriaceae* by the color.

#### 2 EASY READING

The intense colored colonies on a clear agar background facilitate the reading, contrary to VRBG where the smoked background allows little contrast.

### Medium Description

Powder Base	Total ..... 42.3 g/L Agar ..... 11.0 Peptones and Yeast extract ..... 22.0 Mineral salts ..... 4.9 Chromogenic and selective mix ..... 0.7 Growth factors ..... 3.7 Storage at 15/30 °C - pH: 7.4 +/- 0.2 Shelf Life ..... 2 years
Optional Supplements for ready-to-use bottles (to be ordered separately)	1 <sup>st</sup> : Liquid form ..... 1 mL/L Storage at 2-8 °C ..... Shelf Life ..... 2 years 2 <sup>nd</sup> : Powder form qsf 500 mL ..... 0.01 g/L Storage at 2-8 °C ..... Shelf Life ..... 2 years

Usual Samples	Food, foodstuff, environmental samples
Procedure	Pouring, streaking or spreading techniques. Incubation 24 h, 30 °C or 37 °C. Aerobic conditions.

### Order References

Please use these product references when contacting your local distributor:

5000 ml pack ..... EB042

Supplement 1 (to be used only with ready-to-use bottles)..... TT100L

Supplement 2 (to be used only with ready-to-use bottles)..... CV702

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