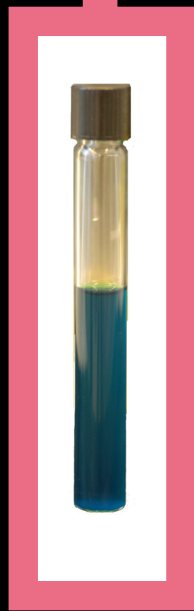
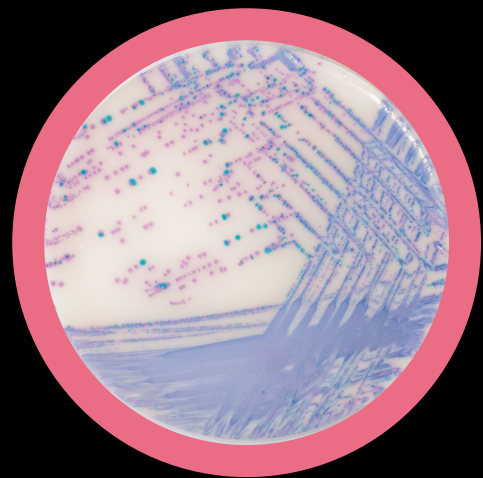
RambaQUICK™ Salmonella Method

Broth enrichment
RambaQUICK™ Salmonella

+

CHROMagar™ Salmonella Plus
Culture media

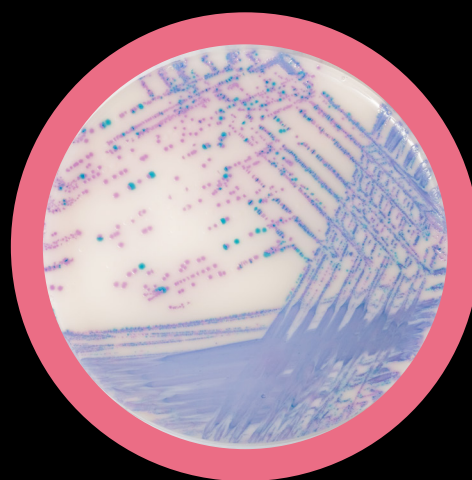
Click below:

● RambaQUICK™ Salmonella Method


Broth enrichment
RambaQUICK™ Salmonella

CHROMagar™ Salmonella Plus
Culture media

For Rapid detection of *Salmonella* species
including *S. typhi*, *S. paratyphi* and
lactose positive *Salmonella*

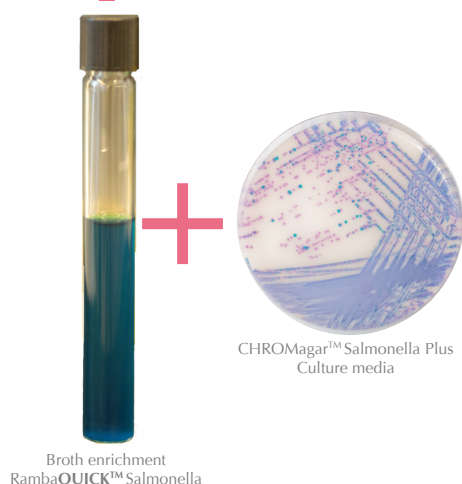www.CHROMagar.com

For Rapid detection of *Salmonella* species including *S. typhi*, *S. paratyphi* and lactose positive strains in food samples

Background

From the CDC report Trends in Foodborne Illness, 1996-2010:

« This year's report (released in CDC's Morbidity and Mortality Weekly Report [MMWR]) summarizes 2010 data from FoodNet and provides the best measure of trends in foodborne disease from 1996–2010. It confirms that *Salmonella* infection has not declined in 15 years,... » One of the report recommendations to reduce the number of *Salmonella* infections is « Enhancing laboratory testing and disease reporting to more quickly identify outbreaks and their causes »

RambaQUICK™ Salmonella Method: Rapid and Simple !

The RambaQUICK™ Salmonella method was designed with the aim of **simplifying the test procedure while maintaining a high level of sensitivity within a conventional culture methodology.**

① TWO IN ONE ENRICHMENT STEP

The Ramba**QUICK**[™] selective broth is a combination of the best functions found in each of both conventional enrichments, the RVS and the MKTTn broths. After the revivification phase in BPW, the *Salmonella* replicates at the exponential growth phase in the optimised Ramba**QUICK**[™] *Salmonella* broth, which offers not only a highly nutrient environment but also inhibits the growth of competitive flora.

② FAST METHOD

- negative in 2 days.
- **positive** in 3 days.

③ SIMPLE

- one selective enrichment.
- one culture plate.

4 MEETS ISO 6579 REQUIREMENTS

by detecting lactose positive *Salmonella* in intense colour on CHROMagar™ *Salmonella* Plus.

- *Salmonella* including *S. typhi*, *S. paratyphi* A and lactose positive *Salmonella* → Mauve

Method Description


RambaQUICK™ Enrichment Broth Powder	Total 34.2 g/L Salts 8.6 Peptone 4.5 Selective mix 21.1 Powder Storage at 15/30°C - pH: 5.7 +/-0.2 Shelf Life: 2 years	
 CHROMagar™ Salmonella Plus Culture Media Powder	Powder Base : Total 32.8 g/L Agar 15.0 Peptone and Yeast extract ... 8.0 Salts 8.5 Chromogenic mix 1.3 Storage at 15/30 °C - pH: 7.5 ± 0.2 Shelf Life: 3 years	Supplements 1st : Liquid form 6.0 Storage at 15/30 °C Shelf Life: 5 years <u>In order to obtain a white opaque background:</u> 2nd : White Opaque Supplement (Powder): 1.0 Storage at 15/30 °C Shelf Life: 3 years

Scientific Publications on this product: available on www.CHROMagar.com

Please read carefully the instructions for use (IFU document) available on www.CHROMagar.com

RambaQUICK™ Salmonella Method

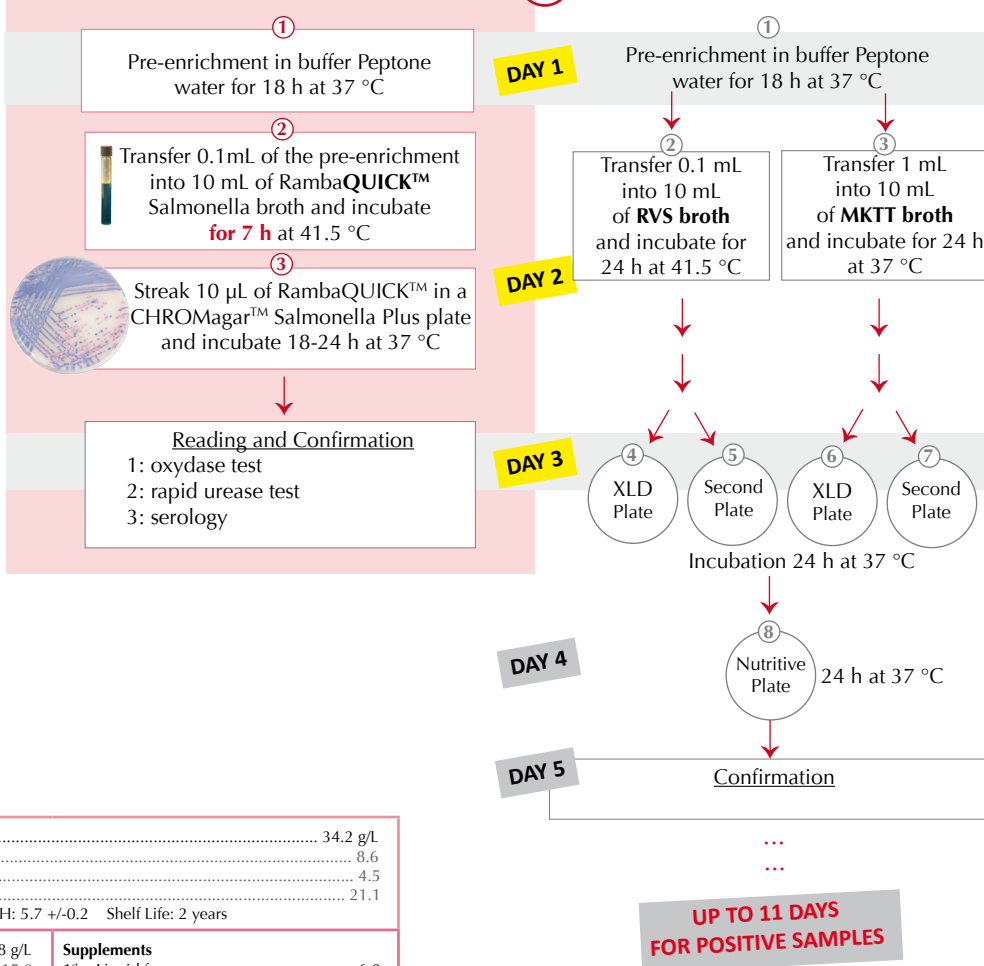VS ISO standard method ISO 6579

**UP TO 11 DAYS
FOR POSITIVE SAMPLES**

F-EXT-036-EN	Version 4.0 / Ago - 25
--------------	------------------------

Order References

Please use these product references when contacting your local distributor:

Ramba**QUICK**TM Salmonella Enrichment broth powder in 1 000 mL pack (100 tests of 10 mL): SQ001
CHROMagarTM Salmonella Plus culture media powder in 5 000 mL pack (250 plates of 20 mL): SA162

RambaQUICK™ Salmonella Method

to prepare 5 000 mL of media (250 tests): 3 packs of SQ001 + 1 pack of SA162

Manufacturer: CHROMagar, 29 avenue George
Sand, 93210 La Plaine Saint-Denis - France

Email: CHROMagar@CHROMagar.com

Website: www.CHROMagar.com

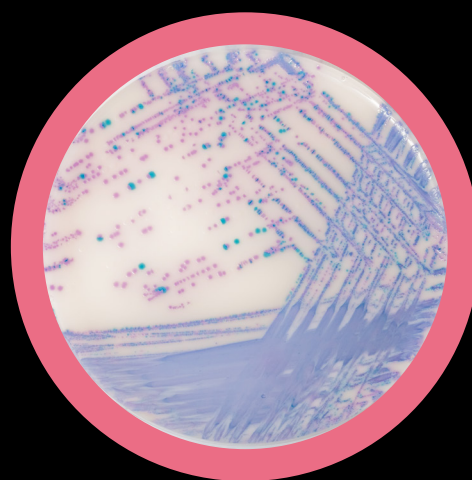
Find your nearest distributor on

Find your nearest distributor on
www.CHROMagar.com/contact

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