# RambaQUICK™ Salmonella

## Instructions For Use

Available in several languages

NT-EXT-025

Version 6.0

Click below for:

ENGLISH

FRANCAIS

ESPAÑOL

DEUTSCH

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RambaQUICK™ Salmonella Broth

RambaQUICK™ Salmonella  
is sold in powder form

# RambaQUICK™ Salmonella

Instructions For Use

ENGLISH

Instructions For Use

TO BE USED WITH CHROMagar™ Salmonella Plus

## MEDIUM PURPOSE

Enrichment broth for rapid detection of *Salmonella* species.

From the CDC report Trends in Foodborne Illness, 1996-2010: « This year's report (released in CDC's Morbidity and Mortality Weekly Report [MMWR]) summarizes 2010 data from FoodNet and provides the best measure of trends in foodborne disease from 1996–2010. It confirms that *Salmonella* infection has not declined in 15 years,... »

One of the report recommendations to reduce the number of *Salmonella* infections is "Enhancing laboratory testing and disease reporting to more quickly identify outbreaks and their causes ».

## COMPOSITION

The product is composed of a powder base

| Product         | = | Base (B)                                       |
|-----------------|---|--|
| Total g/L       |   | 34.2 g/L                                       |
| Composition g/L |   | Peptone 4.5<br>Salts 8.6<br>Selective mix 21.1 |
| Aspect          |   | Powder Form                                    |
| STORAGE         |   | 15-30 °C                                       |
| FINAL MEDIA pH  |   | 5.6 +/- 0.2                                    |

## PREPARATION (Calculation for 1 L)

### Step 1

Preparation of the tubes

- Disperse slowly 34.2 g of powder base in 1 L of purified water.

- Swirl or stir gently to homogenize.

- Heat and bring to boil (100 °C) for 1-2 min while swirling or stirring regularly.

DO NOT HEAT TO MORE THAN 100 °C.

**Warning 1:** If using an autoclave, do so without pressure.

- Pour 10 mL into sterile tubes.

**Advice:** Prepared tubes can be kept for 6 months under refrigeration (2/8 °C) if properly prepared and protected from light.

**Warning 2:** Some crystals could appear later on but do not impact the performances.

## INOCULATION


Remove the tubes from the refrigerator and allow to warm to room temperature before inoculation.

**Please follow the method below.**


### Typical Samples

e.g. food, animal stuff, primary production stage, environmental samples


## METHOD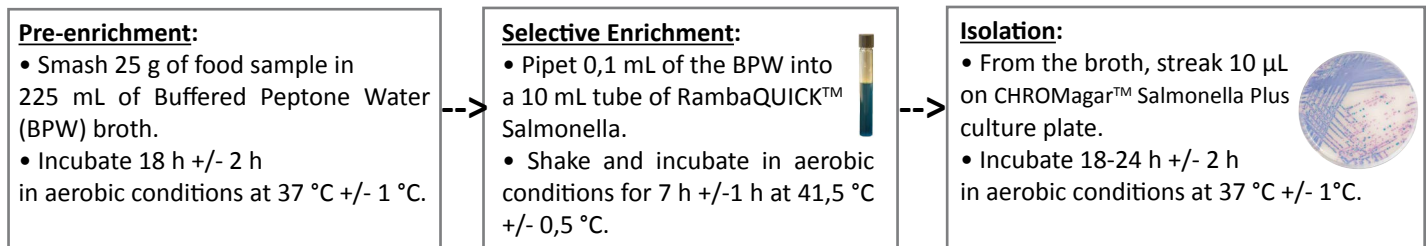

**Typical** colony appearance on CHROMagar™ Salmonella Plus

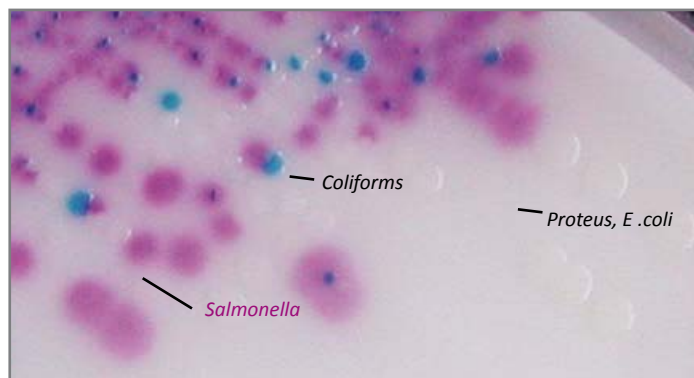The pictures shown are not contractual.
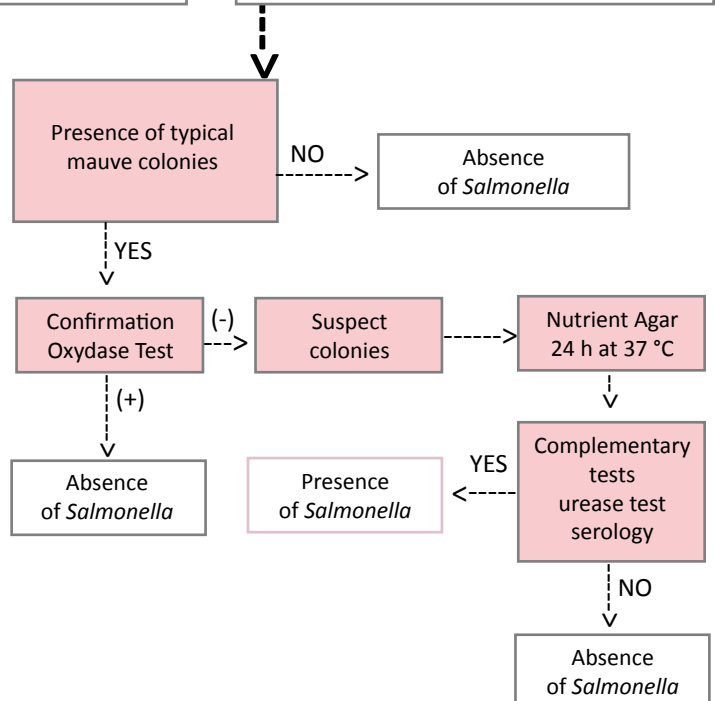
## PERFORMANCE & LIMITATIONS

• Performances of this test depend on chosen *Salmonella* culture medium. We suggest the use of the CHROMagar™ Salmonella Plus medium which allows detection of all *Salmonella* species including *Salmonella typhi*, *S. paratyphi* and lactose positive *Salmonella* according to ISO norm 6579.

## WARNINGS

- Do not use tubes if they show any evidence of contamination or any sign of deterioration.
- Do not use the product beyond its expiry date or if product shows any evidence of contamination or any sign of deterioration.
- For laboratory use. This laboratory product should be used only by trained personnel in compliance with good laboratory practices.
- Any change or modification in the procedure may affect the results.
- Any change or modification of the required storage temperature may affect the performance of the product.
- Unappropriate storage may affect the shelf life of the product.
- Recap the bottles tightly after each preparation and keep them in a low humidity environment, protected from moisture and light.
- For a good microbial detection: collection and transport of specimen should be well handled and adapted to the particular specimen according to good laboratory practices.

## REFERENCES

Please refer to our website page «Publications» for scientific publications about this particular product.

Web link: <http://www.chromagar.com/publication.php>

## DISPOSAL OF WASTE

After use, all tubes and any other contaminated materials must be sterilized or disposed of by appropriate internal procedures and in accordance with local legislations. tubes can be destroyed by autoclaving at 121 °C for at least 20 minutes.

## IFU/LABEL INDEX

-  Catalogue reference
-  Consult instructions for use
-  Quantity of powder sufficient for X liters of media
-  Expiry date
-  Required storage temperature
-  Store away from humidity
-  Protect from light
-  Manufacturer

 Pack Size

1000 mL

100 broths  
of 10 mL

=

Ordering References

SQ001

Weight: 34,2 g

### Need some Technical Documents?

Available  
for download on  
[www.CHROMagar.com](http://www.CHROMagar.com)

- Certificate of Analysis (CoA) --> One per Lot
- Material Safety Data Sheet (MSDS)

CHROMagar™ and Rambach™ are trademarks created by Dr A. Rambach  
ATCC® is a registered trademark of the American Type Culture Collection  
[NT-EXT-025 V6.0 / EN 14-Dec-18](#)

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